

Mouse model for brain metastasis


2019.11.15 Mi Reu Jeong

Materials


1. Xylazine-ketamine for anesthesia
2. Surgical instruments
3. Diameter: 0.38mm polyethylene tubing for catheter
4. 25 μ L Hamilton syringe
5. 4-0 black silk
6. 4-0 nylon
7. Absorbent swabs
8. Blood vessel clamp
9. 10mL syringe
10. Insulin syringe
11. Alcohol
12. PBS
13. 1×10^5 LLC-LUC cells/10 μ L
14. Syringe pump

Methods

1. Connect the catheter to 25 μ L Hamilton syringe
2. Put PBS in a 10mL syringe

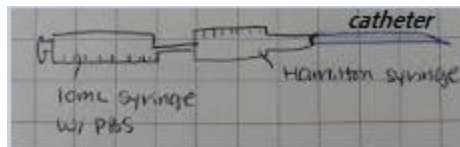

3. Wash 25 μ L Hamilton syringe with PBS
4. Draw 2.5 μ L of air bubble with 25 μ L Hamilton syringe
5. Draw 12.5 μ L of PBS with the same 25 μ L Hamilton syringe
6. Anesthetized a mouse with xylazine-ketamine
7. Expose common carotid artery

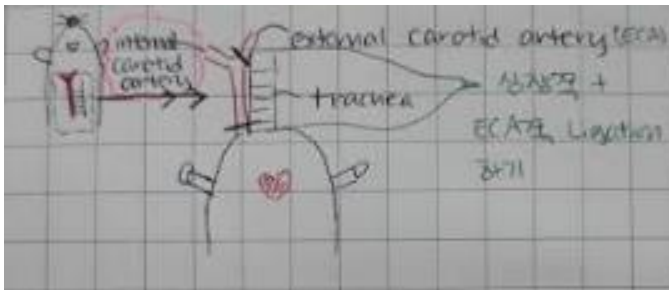


8. Ligate very lower part of common carotid artery
9. Ligate the external carotid artery (the one that looks like directing trachea)



10. Clamp the internal carotid artery



11. Draw $10\mu\text{L}$ of 1×10^5 LLC-LUC cells with $25\mu\text{L}$ the same Hamilton syringe
12. Make a hole on the common carotid artery with insulin syringe
13. Insert the catheter into the lumen of the common carotid artery
14. Tie around the catheter-inserted common carotid artery with black silk
15. Remove the clamp

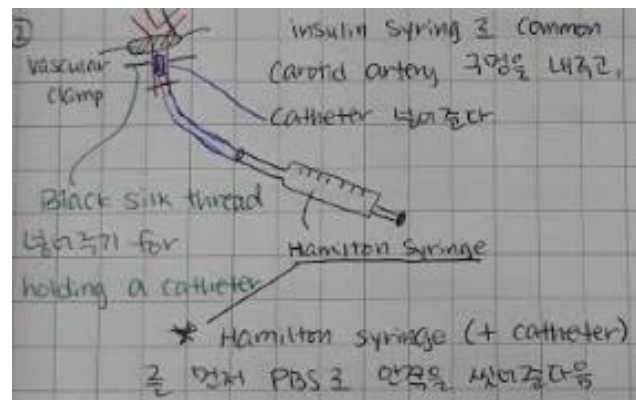

16. Inject the tumor cells 10 μ L/3min. using syringe pump
17. Chase the tumor cells with 12.5 μ L of PBS using syringe pump



18. Remove the catheter
19. Ligate the common carotid artery



20. Suture the opening site with 4-0 nylon
21. Wash the catheter-Hamilton syringe with alcohol

- The brain metastasis would occur at the cerebellum on day 7
- Mouse starts to die on day 10

